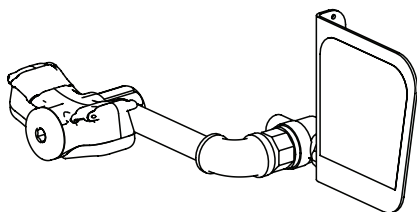
Installation

S19-244 & S19-244FW


Unmounted Halo™ Eyewash & Eye/Face Wash for Laboratory Application

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Model S19-244

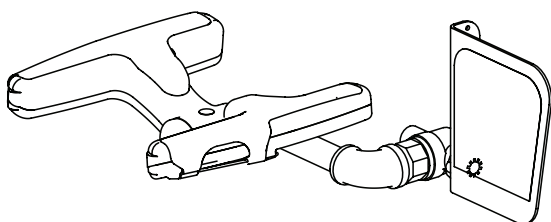Model S19-244FW

Read the instructions in this manual before beginning installation. Save these instructions and refer to them for inspection, maintenance, and troubleshooting information.

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For questions regarding the operation, installation or maintenance of this product, visit bradleycorp.com or call 800.BRADLEY (800.272.3539).

Product warranties and parts information may also be found under "Resources" on our website at bradleycorp.com.

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Table of Contents

Safety Information	3
Supplies Required	3
Assemble Components	4
Connect Water Supply	4
Dimensions	4
Assembly of Components	5
Parts List	5

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To ensure proper operation:

Installation

This fixture has been designed and 3rd party tested to meet ANSI/ISEA Z358.1. The system must be tested periodically to be sure it is functioning properly to meet the standard. Please refer to the ANSI/ISEA Z358.1 standard on testing requirements.

The installation, maintenance and location of all eyewash and eye/face washes must comply with the requirements of ANSI/ISEA Z358.1. Workers who may come in contact with potentially hazardous materials should be trained regarding the location and proper operation of emergency equipment per ANSI/ISEA Z358.1.

ANSI/ISEA Z358.1 requires this unit to be used with a clean, potable, uninterrupted supply of tepid water. Bradley plumbed emergency fixtures require a minimum of 30 psi flowing pressure in order to function as designed and meet the ANSI/ISEA Z358.1 standard.

Before installation make sure that this equipment will meet the requirements of the potential hazardous contaminants in your location.

Installation of this system must be completed by a qualified plumber in compliance with all national and local codes. Compliance and conformity to drain requirements and other local codes and ordinances is the responsibility of the installer.

Weekly Inspections

To ensure all eyewash and eye/face wash fixtures are ready to be used when needed, regular checks and maintenance must be conducted. Weekly inspections must be conducted on all safety equipment to make sure a suitable flushing fluid supply is present and to make sure the equipment is in good operating condition.

Water Temperature

ANSI/ISEA Z358.1 requires tepid water. Suitable range is 60°F to 100°F (16°C to 38°C). Personal injury is possible outside this temperature range.

Supplies Required

- Pipe sealant
- Piping to 1/2" NPT water supply inlet on unit
- Sign-mounting hardware
- Backflow prevention device as required by UPC, IPC and/or local plumbing codes

Local codes may require the installation of a backflow prevention valve to complete proper installation. Backflow prevention valves are not included with the fixture and may be supplied by the contractor or purchased from Bradley Corporation.

Assemble Components

 The top of the eyewash heads should be 33–45" (838–1143 mm) from the floor.

1. Select the desired location for the eyewash on the sink deck and drill a 7/8" diameter hole.
2. Assemble the components as shown on page 5.
 - Apply pipe sealant (supplied by installer) to all male-threaded pipe joints.
 - Use a strap wrench around pipes when tightening to prevent marring the BradTect® protective, safety yellow coating.

Connect Water Supply

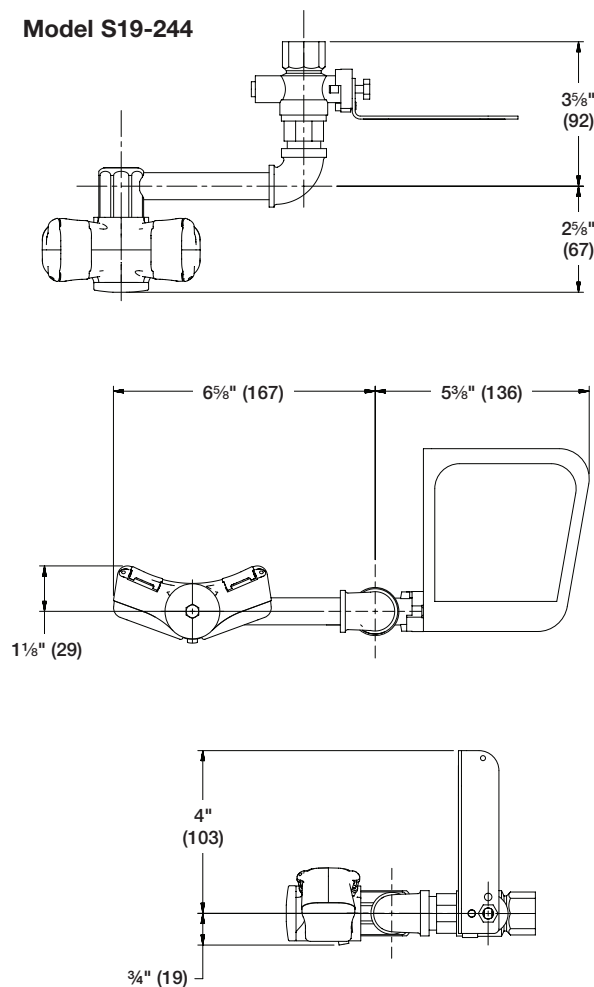
1. Connect the water supply piping (by installer) to the 1/2" NPT inlet on the eyewash.
2. Open the water supply lines. Test for leaks and adequate water flow.
3. Mount the safety sign to the wall using the sign-mounting hardware supplied by the installer.

 All dimensions assume standard thread engagement. Variations in manufacturing allow for $\pm 1/8$ " (3 mm) per thread joint. To find the tolerance of a dimension, add the number of thread joints in between a dimension and multiply it by 1/8" (3 mm). Connect the water supply using installer supplied supply line. Mount the safety sign to wall or pipe using sign-mounting hardware supplied by installer. Open the water supply lines. Test for leaks and adequate water flow.


Dimensions

(mm)


Model S19-244

Model S19-244FL

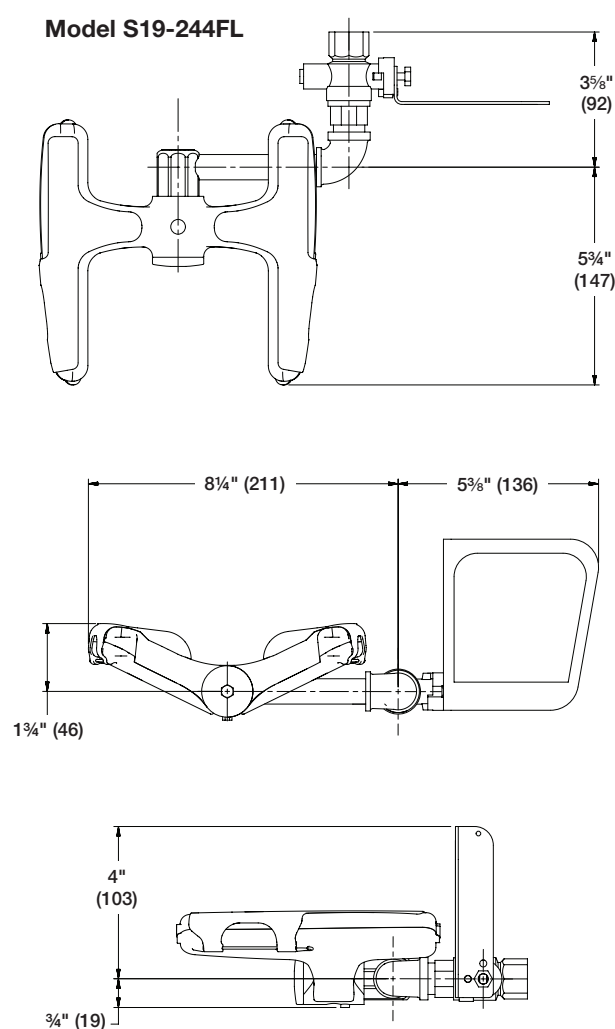Assembly of Components

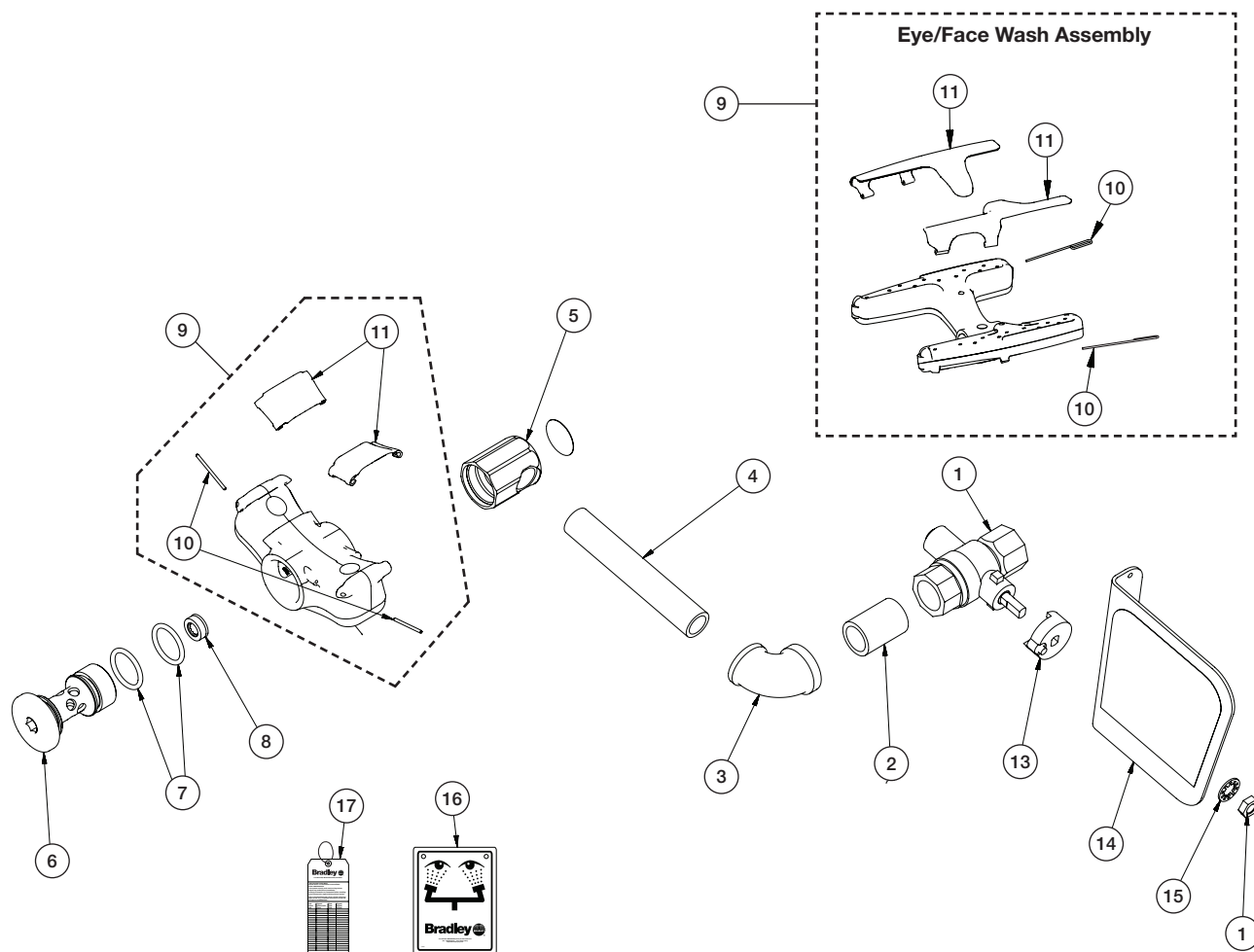Parts List

Item	Part No.	Qty.	Description
1	S27-327	1	1/2" Ball Valve, Brass
2	113-006AW	1	Pipe 1/2" X Close Chrome Brass Red Brass SCH 40
3	169-1023A	1	1/2" X 3/8" Red Elbow Chrome
4	113-1192	1	Pipe 3/8" X 4" Chrome Red Brass SCH 40
5	171-005	1	Body Mount - Short 3/8 NPT
6	115-169	1	Body Cap - Short
7	125-157	2	O-Ring (-117), .799 ID X .103 W EPDM- 70 DURO-EU-CERT
8	125-175	1	Halo Flow Control
9	S05-190	1	Sprayhead Assembly - Eyewash (Includes Items 10 & 11) (S19-244)
	S05-192	1	Sprayhead Assembly - Eye/Face Wash (Includes Items 10 & 11) (S19-244FW)

Item	Part No.	Qty.	Description
10	Included in S05-190	2	Hinge Clip (S19-244)
	Included in S05-192	2	Hinge Clip (S19-244FW)
11	Included in S05-190	2	Dust Cover - Eyewash (S19-244)
	Included in S05-192	1	Dust Cover - Eye/Face Wash - Right (S19-244FW)
	Included in S05-192	1	Dust Cover - Eye/Face Wash - Left (S19-244FW)
13	153-372R	1	Ball Valve Adaptor - RH
14	S08-336	1	RH Handle / Label Assembly
15	142-002DA	1	Washer 5/16 Int-Tooth Lock 18-8-SS
16	114-051	1	Safety Sign Eyewash
17	204-421	1	Emergency Inspection Tag

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© 2023 Bradley